



vista[®]
Patent Pending

for...

**LIQUID DETERGENT
(CLEAR)**



Major brand owners know the importance that color appeal and color consistency play in a consumer's decision to purchase their Liquid Detergent products. Vista ensures that the color and clarity of Liquid Detergents conform to quality standards by simultaneously measuring both transmittance color and haze% in a single measurement. Any issues can be immediately isolated, the root cause discovered, and the process corrected.

Vista includes easy to navigate embedded EasyMatch Essentials quality control software and modern data streaming and communications, including USB, Ethernet and Wi-Fi. With its capacitive high-resolution color touchscreen display, Vista sets a new standard for transmission color and appearance analysis of water products.



**Simultaneously Measure
Transmission Color and Haze%
in a Single Measurement**



**Long-life, Full Spectrum LED
Illumination Guaranteed for
Five Years**



**One-touch standardization -
no standards or accessories
required**



**Embedded EasyMatch
Essentials Quality Control
Software – No PC Required**



**Smart User Interface and Data
Management – USB, Ethernet
and WiFi**




measure
color
the way your
eye sees it

The Best Solution for Color Measurement of Clear Liquid Detergent




Simultaneously Measure Transmittance Color and Haze%

Simultaneously measure transmittance color and haze% compliant with ASTM D1003 Procedure B, in a single measurement, along with a host of other color scales and color indices - in under seven seconds.

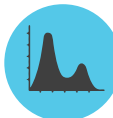


Full Spectrum, Balanced LED Illumination

Illuminating from 400-700 nm, the balanced, full spectrum LED illumination is guaranteed to perform to original factory specifications for a minimum of 5 years of continuous measurements for years of worry-free operation.

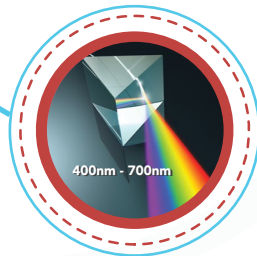

Operator-friendly One-touch Standardization

Instrument standardization is fully automated, no tiles or accessories are required. Simply press the 'standardize' icon, and the instrument automatically standardizes for both color and haze.

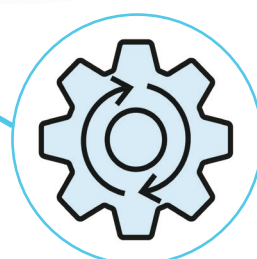

Embedded EasyMatch Essentials Quality Control Software – No PC Required

Simultaneously Measure transmittance color and Haze% compliant with ASTM D1003 Procedure B, in a single measurement, along with a host of other color scales and color indices in Vista's embedded EasyMatch Essentials quality control software.



Smart User Interface and Data Management

View measurement results and color graphics on the 7-inch 1280x800 capacitive industrial touchscreen, and easily communicate data where it needs to go via USB, USB on-the-go, Ethernet, Wi-Fi, and built in email.

EasyMatch Essentials

Keys for the Successful Measurement of Clear Liquid Detergent



D02-1017-192
Multi-Function Sample Holder



D02-1017-344
Macro Cell Holder for Color
and Haze Measurement

*HunterLab provides a variety of standard cell holder options.
Contact us for help in selecting the best solution to fit your needs.*

L02-1018-683
10 mm Flow Through Cell Holder Kit
Includes:

- 10 mm Flow Through Cell (C04-1001-958)
- Cell Holder Assembly (D02-1018-678)
- 1 meter, 7 mm ID tubing
- (4) Flow Control Chokes (Various apertures) (L02-1018-101)

20 mm Flow Through Cell Kit also available

